

# *PRODUCT INFORMATION AND MANUAL*

## ***mouse TNF- $\alpha$ Instant ELISA***

### ***BMS607INST***

Enzyme-linked immunosorbent assay for quantitative  
detection of mouse TNF- $\alpha$ .

For research use only.

Not for diagnostic or therapeutic procedures.

▽ 128 Tests



**mouse TNF- $\alpha$   
BMS607INST**



Bender MedSystems GmbH  
Campus Vienna Biocenter 2  
A-1030 Vienna, Austria, Europe



## TABLE OF CONTENTS

1	Intended Use	4
2	Summary	4
3	Principles of the Test	5
4	Reagents Provided	6
5	Storage Instructions	6
6	Specimen Collection	7
7	Materials Required But Not Provided	8
8	Precautions for Use	9
9	Preparation of Reagents and Samples	11
10	Test Protocol	12
11	Calculation of Results	16
12	Limitations	19
13	Performance Characteristics	20
14	Bibliography	22
15	Ordering Information	23
16	Reagent Preparation Summary	24
17	Test Protocol Summary	25

## 1 Intended Use

The mouse TNF- $\alpha$  Instant ELISA is an enzyme-linked immunosorbent assay for the quantitative detection of mouse TNF- $\alpha$ . The mouse TNF- $\alpha$  Instant ELISA is **for research use only. Not for diagnostic or therapeutic procedures.**

## 2 Summary

Tumor Necrosis Factor  $\alpha$  (TNF $\alpha$ ), also known as cachectin, is a polypeptide cytokine produced by monocytes and macrophages. It functions as a multipotent modulator of immune response and further acts as a potent pyrogen (1, 4). TNF $\alpha$  circulates throughout the body responding to stimuli (infectious agents or tissue injury), activating neutrophils, altering the properties of vascular endothelial cells, regulating metabolic activities of other tissues, as well as exhibiting tumoricidal activity by inducing localized blood clotting. TNF $\alpha$  also inhibits lipoprotein lipase activity resulting in cachexia, a physical wasting condition (1, 4). Due to its varied actions throughout the immune system, TNF $\alpha$  may play a role in the pathogenesis of many disease states.

TNF $\alpha$  production is mediated by the action of lymphokines and endotoxins on the macrophage. Purified monocytes produce TNF $\alpha$  within four hours of stimulation by recombinant IL-2 (3) and there is some *in vitro* evidence to suggest that TNF $\alpha$  is expressed at high levels and with prolonged kinetics in T cells stimulated by both CD2 and CD28 (2). Secretion of TNF $\alpha$  is enhanced by gamma interferon. TNF then induces or enhances the specific production of Class I MHC antigen, GM-CSF, and IL-1. Recent evidence has suggested an intracellular role for this peptide (5).

### 3 Principles of the Test

An anti-mouse TNF- $\alpha$  coating antibody is adsorbed onto microwells. Mouse TNF- $\alpha$  present in the sample or standard binds to antibodies adsorbed to the microwells; a biotin-conjugated anti-mouse TNF- $\alpha$  antibody binds to mouse TNF- $\alpha$  captured by the first antibody. Streptavidin-HRP binds to the biotin conjugated anti-mouse TNF- $\alpha$ .

Following incubation unbound biotin conjugated anti-mouse TNF- $\alpha$  and Streptavidin-HRP is removed during a wash step, and substrate solution reactive with HRP is added to the wells.

A coloured product is formed in proportion to the amount of soluble mouse TNF- $\alpha$  present in the sample. The reaction is terminated by addition of acid and absorbance is measured at 450 nm. A standard curve is prepared from 7 mouse TNF- $\alpha$  standard dilutions and mouse TNF- $\alpha$  sample concentration determined.

Figure 1

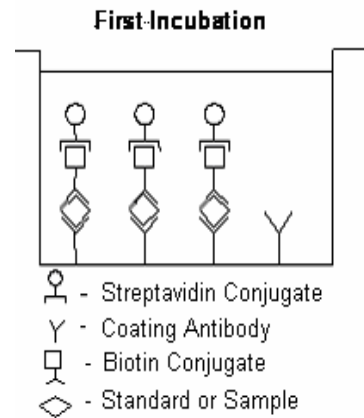


Figure 2


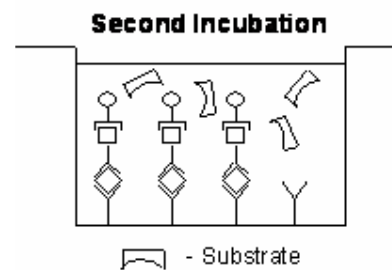


Figure 3

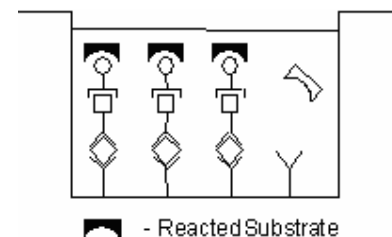

#### 4 Reagents Provided

- 1 aluminium pouch with a **Microwell Plate coated with polyclonal Antibody** to mouse TNF- $\alpha$ , **Biotin-Conjugate** (anti-mouse TNF- $\alpha$  polyclonal antibody), **Streptavidin-HRP** and Sample Diluent, lyophilized
- 2 aluminium pouches with a mouse TNF- $\alpha$  **Standard curve** (coloured)
- 1 bottle (25 ml) **Wash Buffer Concentrate** 20x (phosphate-buffered saline with 1% Tween 20)
- 1 vial (12 ml) **Sample Diluent** (Use when an external predilution of the samples is needed)
- 1 vial (15 ml) **Substrate Solution** (tetramethyl-benzidine)
- 1 vial (15 ml) **Stop Solution** (1M Phosphoric acid)
- 2 **Adhesive Films**

#### 5 Storage Instructions

Store ELISA plate and Standard curves or whole kit at -20°C. The plate and the standard can also be removed, stored at -20°C, remaining kit reagents can be stored between 2° and 8°C. Expiry of the kit and reagents is stated on labels.

The expiry of the kit components can only be guaranteed if the components are stored properly, and if, in case of repeated use of one component, the reagent is not contaminated by the first handling.

## 6 Specimen Collection

Serum and cell culture supernatant samples are suitable for use in the assay. Remove serum from the clot as soon as possible after clotting.

Samples containing a visible precipitate must be clarified prior to use in the assay. Do not use grossly hemolyzed or lipemic specimens.

Samples must be stored frozen at  $-20^{\circ}\text{C}$  to avoid loss of bioactive mouse TNF- $\alpha$ . If samples are to be run within 24 hours, they may be stored at  $2^{\circ}$  to  $8^{\circ}\text{C}$  (for sample stability refer to 13).

Avoid repeated freeze-thaw cycles. Prior to assay, frozen serum or plasma should be brought to room temperature slowly and mixed gently.

## 7 Materials Required But Not Provided

- 5 ml and 10 ml graduated pipettes
- 5  $\mu$ l to 1000  $\mu$ l adjustable single channel micropipettes with disposable tips
- 50  $\mu$ l to 300  $\mu$ l adjustable multichannel micropipette with disposable tips
- Multichannel micropipette reservoir
- Beakers, flasks, cylinders necessary for preparation of reagents
- Device for delivery of wash solution (multichannel wash bottle or automatic wash system)
- Microwell strip reader capable of reading at 450 nm (620 nm as optional reference wave length)
- Glass-distilled or deionized water
- Statistical calculator with program to perform linear regression analysis

## 8 Precautions for Use

- All chemicals should be considered as potentially hazardous. We therefore recommend that this product is handled only by those persons who have been trained in laboratory techniques and that it is used in accordance with the principles of good laboratory practice. Wear suitable protective clothing such as laboratory overalls, safety glasses and gloves. Care should be taken to avoid contact with skin or eyes. In the case of contact with skin or eyes wash immediately with water. See material safety data sheet(s) and/or safety statements(s) for specific advice.
- Reagents are intended for research use only and are not for use in diagnostic or therapeutic procedures.
- Do not mix or substitute reagents with those from other lots or other sources.
- Do not use kit reagents beyond expiration date on label.
- Do not expose kit reagents to strong light during storage or incubation.
- Do not pipette by mouth.
- Do not eat or smoke in areas where kit reagents or samples are handled.
- Avoid contact of skin or mucous membranes with kit reagents or specimens.
- Rubber or disposable latex gloves should be worn while handling kit reagents or specimens.
- Avoid contact of substrate solution with oxidizing agents and metal.
- Avoid splashing or generation of aerosols.
- In order to avoid microbial contamination or cross-contamination of reagents or specimens which may invalidate the test use disposable pipette tips and/or pipettes.
- Use clean, dedicated reagent trays for dispensing substrate reagent.

- Glass-distilled water or deionized water must be used for reagent preparation.
- Substrate solution must be at room temperature prior to use.
- Decontaminate and dispose specimens and all potentially contaminated materials as they could contain infectious agents. The preferred method of decontamination is autoclaving for a minimum of 1 hour at 121.5°C.
- Liquid wastes not containing acid and neutralized waste may be mixed with sodium hypochlorite in volumes such that the final mixture contains 1.0% sodium hypochlorite. Allow 30 minutes for effective decontamination. Liquid waste containing acid must be neutralized prior to the addition of sodium hypochlorite.

## 9 Preparation of Reagents and Samples

The buffer concentrate should be brought to room temperature and diluted before starting the test procedure.

### 9.1 Wash Buffer

If crystals have formed in the **Wash Buffer Concentrate**, warm it gently until they have completely dissolved.

Pour entire contents (25 ml) of the Wash Buffer Concentrate into a clean 500 ml graduated cylinder. Bring to final volume to 500 ml with glass-distilled or deionized water. Mix gently to avoid foaming. The pH of the final solution should adjust to 7.4.

Transfer to a clean wash bottle and store at 2° to 25°C. Please note that Wash Buffer is stable for 30 days.

## 10 Test Protocol

- **Use plate immediately after removal from -20°C!**
  - **Do not wait until pellets have completely dissolved before applying samples - the binding reaction in the standard strips starts immediately after addition of water!**
  - **Do not try to dissolve pellets by pipetting up and down in the wells - some parts of the pellet could stick to the tip creating high variation of results**
  - **Perform the washing step with at least 400 µl of washing buffer as stated in the manual or fill the wells completely - otherwise any pellet residues sticking to the rim of the well will not be removed and create high variation of results**
  - **Allow the washing buffer to sit in the wells for a few seconds before aspiration**
  - **Remove covers of the standard strips carefully in order that all the lyophilised pellets remain in the wells**
- a. Determine the **number of Microwell Strips** required to test the desired number of samples plus Microwell Strips for blanks and standards (coloured). Each sample, standard and blank should be assayed in duplicate. Remove extra Microwell Strips from holder and store in foil bag with the desiccant provided at -20°C sealed tightly. Place Microwell Strips containing the standard curve in position A1/A2 to H1/H2 (see Table 1).
- b. Add 100 µl of **distilled water** to the sample wells.
- c. Add **distilled water** to all standard and blank wells as indicated on the label of the standard strips (A1, A2 to H1, H2).

Table 1

Table depicting an example of the arrangement of blanks, standards and samples in the microwell strips:

	<b>1</b>	<b>2</b>	<b>3</b>	<b>4</b>
<b>A</b>	Standard 1 (2000 pg/ml)	Standard 1 (2000 pg/ml)	Sample 1	Sample 1
<b>B</b>	Standard 2 (1000 pg/ml)	Standard 2 (1000 pg/ml)	Sample 2	Sample 2
<b>C</b>	Standard 3 (500 pg/ml)	Standard 3 (500 pg/ml)	Sample 3	Sample 3
<b>D</b>	Standard 4 (250 pg/ml)	Standard 4 (250 pg/ml)	Sample 4	Sample 4
<b>E</b>	Standard 5 (125 pg/ml)	Standard 5 (125 pg/ml)	Sample 5	Sample 5
<b>F</b>	Standard 6 (62,5 pg/ml)	Standard 6 (62,5 pg/ml)	Sample 6	Sample 6
<b>G</b>	Standard 7 (31,3 pg/ml)	Standard 7 (31,3 pg/ml)	Sample 7	Sample 7
<b>H</b>	Blank	Blank	Sample 8	Sample 8

- d. Add 50  $\mu$ l of each **Sample**, in duplicate, to the designated wells and mix the contents.
- e. Cover with a **adhesive film** and incubate at room temperature (18°C to 25°C) for 3 hours, if available on a microplate shaker at 200 rpm.
- f. Remove adhesive film and empty wells. **Wash** the microwell strips 6 times with approximately **400  $\mu$ l** Wash Buffer per well with thorough aspiration of microwell contents between washes. Allow the Wash Buffer to sit in the wells for about **10 – 15 seconds** before aspiration. Take care not to scratch the surface of the microwells.

After the last wash, tap microwell strips on absorbent pad or paper towel to remove excess Wash Buffer. Use the microwell strips immediately after washing or place upside down on a wet absorbent paper for no longer than 15 minutes. Do not allow wells to dry.

- g. Pipette 100  $\mu$ l of **TMB Substrate Solution** to all wells, including the blank wells.

**The OD values of the standard curve may vary according to the conditions of assay performance** (e.g. operator, pipetting technique, washing technique or temperature effects). Furthermore shelf life of the kit may affect enzymatic activity and thus colour intensity. Values measured are still valid.

- h. Incubate the microwell strips at room temperature (18° to 25°C) for 10 min. Avoid direct exposure to intense light.
- i. Stop the enzyme reaction by quickly pipetting 100  $\mu$ l of **Stop Solution** into each well, including the blank wells. It is important that the Stop Solution is spread quickly and uniformly throughout the microwells to completely inactivate the enzyme. Results must be read within 5 min after the Stop Solution is added.
- j. Read absorbance of each microwell on a spectro-photometer using 450 nm as the primary wave length (optionally 620 nm as the reference wave length; 610 nm to 650 nm is acceptable). Blank the plate reader according to the manufacturer's instructions by using the blank wells. Determine the absorbance of both the samples and the mouse TNF- $\alpha$  standards.

**Note: In case of incubation without shaking the obtained O.D. values may be lower than indicated below. Nevertheless the results are still valid.**

## 11 Calculation of Results

- Calculate the average absorbance values for each set of duplicate standards and samples. Duplicates should be within 20 per cent of the mean.
- Create a standard curve by plotting the mean absorbance for each standard concentration on the ordinate against the mouse TNF- $\alpha$  concentration on the abscissa. Draw a best fit curve through the points of the graph.
- To determine the concentration of circulating mouse TNF- $\alpha$  for each sample, first find the mean absorbance value on the ordinate and extend a horizontal line to the standard curve. At the point of intersection, extend a vertical line to the abscissa and read the corresponding mouse TNF- $\alpha$  concentration.
- It is suggested that each testing facility establishes a control sample of known mouse TNF- $\alpha$  concentration and runs this additional control with each assay. If the values obtained are not within the expected range of the control, the assay results may be invalid.
- A representative standard curve is shown in Figure 4. This curve cannot be used to derive test results. Every laboratory must prepare a standard curve for each group of microwell strips assayed.
- If the sample is prediluted the value read from the standard curve must be multiplied by the dilution factor. For example, if the sample is prediluted 1:25 the value read from the standard curve must be multiplied by a factor of 25 to give the final result.

Figure 4

Representative standard curve for mouse TNF- $\alpha$  Instant ELISA. Mouse TNF- $\alpha$  was diluted in serial 2-fold steps in Sample Diluent each symbol represents the mean of 3 parallel titrations.

Do not use this standard curve to derive test results. A standard curve must be run for each group of microwell strips assayed.

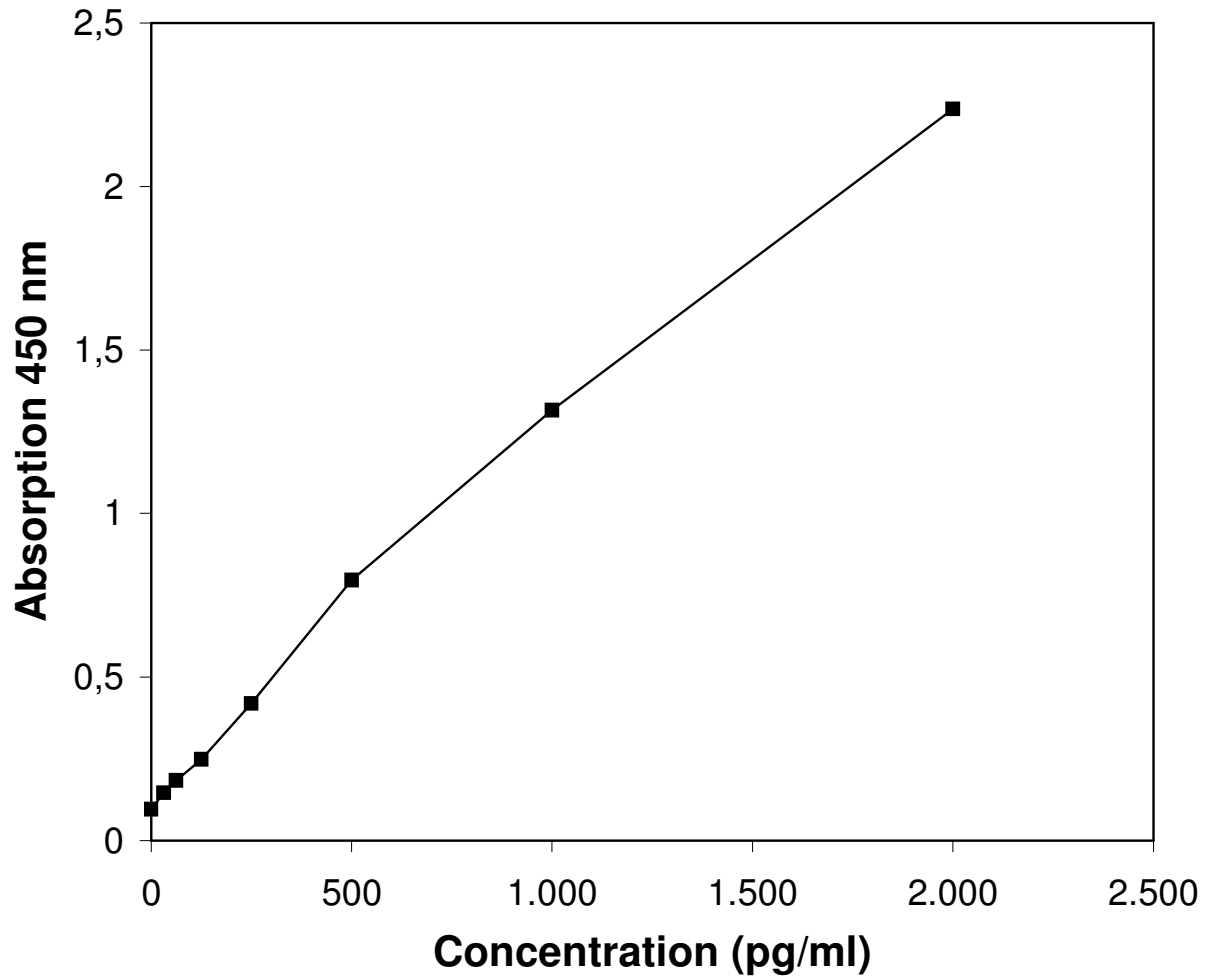


Table 2

Typical data using the mouse TNF- $\alpha$  INSTANT ELISA

Measuring wavelength: 450 nm

Reference wavelength: 620 nm

Standard	mouse TNF $\alpha$ Concentration (pg/ml)	O.D. (450 nm)	O.D. Mean	C.V. (%)
1	2000	2.250	2.237	0.8
	2000	2.224		
2	1000	1.274	1.316	4.5
	1000	1.358		
3	500	0.820	0.796	4.3
	500	0.772		
4	250	0.414	0.419	1.7
	250	0.424		
5	125	0.254	0.249	3.1
	125	0.243		
6	62,5	0.190	0.184	5.0
	62,5	0.177		
7	31,3	0.155	0.147	8.2
	31,3	0.138		
Blank		0.102	0.096	
		0.090		

## 12 Limitations

- Since exact conditions may vary from assay to assay, a standard curve must be established for every run.
- Bacterial or fungal contamination of either screen samples or reagents or cross-contamination between reagents may cause erroneous results.
- Disposable pipette tips, flasks or glassware are preferred, reusable glassware must be washed and thoroughly rinsed of all detergents before use.
- Improper or insufficient washing at any stage of the procedure will result in either false positive or false negative results. Empty wells completely before dispensing fresh wash solution, fill with Wash Buffer as indicated for each wash cycle and do not allow wells to sit uncovered or dry for extended periods.

## **13 Performance Characteristics**

### **13.1 Sensitivity**

The limit of detection of mouse TNF- $\alpha$  defined as the analyte concentration resulting in an absorbance significantly higher than that of the dilution medium (mean plus 2 standard deviations) was determined to be 4,3 pg/ml (mean of 6 independent assays).

### **13.2 Reproducibility**

#### **13.2.1 Intra-assay**

Reproducibility within the assay was evaluated in 3 independent experiments. Each assay was carried out with 6 replicates of 8 serum samples containing different concentrations of mouse TNF- $\alpha$ . 2 standard curves were run on each plate. The calculated overall intra-assay coefficient of variation was < 10%.

#### **13.2.2 Inter-assay**

Assay to assay reproducibility within one laboratory was evaluated in 3 independent experiments. Each assay was carried out with 6 replicates of 8 serum samples containing different concentrations of mouse TNF- $\alpha$ . 2 standard curves were run on each plate. The calculated overall coefficient of variation was < 10%.

### **13.3 Spike Recovery**

The spike recovery was evaluated by spiking 4 levels of mouse TNF- $\alpha$  into serum. Recoveries were determined in 3 independent experiments with 4 replicates each. The unspiked serum was used as blank in these experiments. The overall mean recovery was 98,0 %.

### **13.4 Dilution Parallelism**

4 serum samples with different levels of mouse TNF- $\alpha$  were analysed at serial 2 fold dilutions with 4 replicates each. The overall mean recovery was 105,0 %.

## **13.5 Sample Stability**

### **13.5.1 Freeze-Thaw Stability**

Aliquots of serum samples (unspiked or spiked) were stored at  $-20^{\circ}\text{C}$  and thawed 5 times, and the mouse TNF- $\alpha$  levels determined. There was no significant loss of mouse TNF- $\alpha$  immunoreactivity detected by freezing and thawing.

### **13.5.2 Storage Stability**

Aliquots of serum samples (spiked or unspiked) were stored at  $-20^{\circ}\text{C}$ ,  $2-8^{\circ}\text{C}$ , room temperature (RT) and at  $37^{\circ}\text{C}$ , and the mouse TNF- $\alpha$  level determined after 24h. There was no significant loss of mouse TNF- $\alpha$  immunoreactivity detected during storage at  $-20^{\circ}\text{C}$  and  $4^{\circ}\text{C}$ . Storage at room temperature and  $37^{\circ}\text{C}$  gave rise to loss of mTNF- $\alpha$  immunoreactivity.

## **13.6 Specificity**

The interference of circulating factors of the immune system was evaluated by spiking these proteins at physiologically relevant concentrations into serum. There was no crossreactivity detected.

## **13.7 Hook effect**

Samples with expected concentrations 2 fold higher than the concentration of the highest standard should be diluted 10 fold in Sample Diluent before assay performance to prevent false negative results due to "hook effects".

## **13.8 Expected Values**

A panel of sera samples from randomly selected apparently healthy mice was tested for mouse TNF- $\alpha$ . There were no detectable mouse TNF- $\alpha$  levels found. Elevated mouse TNF- $\alpha$  levels depend on the type of immunological disorder.

## 14 Bibliography

- 1) Beutler B., et al. (1988). *Endo. Rev.* 9, 57.
- 2) Cerdan C., et al. (1991). *J. Immunol.* 146, 560.
- 3) Economou J. S., et al. (1989). *Immunol.* 67, 514.
- 4) Maury C. P. J. (1986). *Acta Med. Scan.* 220, 387.
- 5) Smith M. R., et al. (1990). *J. Immunol.* 144, 560.

For literature update refer to **[www.bendermedsystems.com](http://www.bendermedsystems.com)**

## 15 Ordering Information

For orders please contact:

### **Europe-Headquarters**

Bender MedSystems GmbH  
Campus Vienna Biocenter 2  
A-1030 Vienna, Austria, Europe  
phone: +43 1 796 40 40 ext. 114  
fax: +43 1 796 40 40 ext. 400  
e-mail: [order@bendermedsystems.com](mailto:order@bendermedsystems.com)

### **USA**

Bender MedSystems, Inc.  
849 Hinckley Road  
Burlingame, CA 94010, USA  
toll-free phone: +1 (866) 952 2112  
toll-free fax: +1 (877) 952 2112  
phone: +1 (650) 952 2112  
fax: +1 (650) 952 2252

For technical information please contact:

e-mail: [techserv@bendermedsystems.com](mailto:techserv@bendermedsystems.com)  
**[www.bendermedsystems.com](http://www.bendermedsystems.com)**

**Cat.No. BMS607INST mouse TNF- $\alpha$  INSTANT ELISA**

## **16 Reagent Preparation Summary**

### **16.1 Wash Buffer**

Add **Wash Buffer Concentrate** 20 x (25 ml) to 475 ml distilled water

## 17 Test Protocol Summary

- Place standard strips in position A1/A2 to H1/H2.
- Add 100 µl **distilled water** to sample wells.
- Add **distilled water**, in duplicate, to all standard and blank wells as indicated on the label of the standard strips.
- Add 50 µl **Sample** to designated wells.
- Cover microwell strips and incubate 3 hours at room temperature (18° to 25°C) if available on a microplate shaker at 200 rpm.
- Empty and wash microwell strips 6 times with 400 µl **Wash Buffer**.
- Add 100 µl of **TMB Substrate Solution** to all wells including blank wells.
- Incubate the microwell strips for 10 minutes at room temperature (18° to 25°C).
- Add 100 µl **Stop Solution** to all wells including blank wells.
- Blank microwell reader and measure colour intensity at 450 nm within 5 minutes.